



SA201(-EC) Spectrum Analyzer Controller

Operating Manual



















Table of Contents

Chapter 1	Warning Symbol Definitions	1
Chapter 2	Safety	2
Chapter 3	Overview.....	3
	3.1. Parts List.....	3
	3.2. Compatible Fabry-Perot Scanning Heads	4
Chapter 4	Description.....	5
	4.1. Detector Gain Adjustment.....	6
	4.2. DC Offset Control	6
	4.3. Sweep Expansion Control	6
	4.4. Waveform Control	6
	4.5. Power Switch.....	6
	4.6. Power On Indicator	6
	4.7. Amplitude Control.....	7
	4.8. Rise Time Control	7
	4.9. Trigger Output BNC.....	8
	4.10. Output BNC.....	8
	4.11. Ground Plug.....	9
	4.12. AC Input Connector	9
	4.13. PD Amplifier Input BNC	9
	4.14. PD Amplifier Output BNC.....	10
	4.15. Voltage Selector Switch.....	10
Chapter 5	Operation.....	11
	5.1. PD Blanking Circuit.....	11
	5.2. Replacing the Fuse	12
	5.2.1. Materials Needed.....	12
	5.2.2. Fuse Replacement.....	12

5.3.	<i>Selecting the Line Voltage</i>	13
5.4.	<i>Cleaning</i>	13
Chapter 6	Recommended Setup	14
6.1.	<i>Recommended Setup for SA200 Series Fabry-Perot Interferometers (Except SA200-30C)</i>	14
6.2.	<i>Recommended Setup for SA200-30C Fabry-Perot Interferometer</i>	15
Chapter 7	Specifications	17
7.1.	<i>Output Characteristics</i>	17
7.2.	<i>Trigger Characteristics</i>	17
7.3.	<i>Photodiode Amplifier Characteristics</i>	18
7.4.	<i>Physical Features</i>	19
7.5.	<i>Power Supply</i>	19
Chapter 8	Regulatory	20
Chapter 9	Declaration of Conformity	21
Chapter 10	Thorlabs Worldwide Contacts	22

Chapter 1 Warning Symbol Definitions

Below is a list of warning symbols you may encounter in this manual or on your device.

Symbol	Description
	Direct Current
	Alternating Current
	Both Direct and Alternating Current
	Earth Ground Terminal
	Protective Conductor Terminal
	Frame or Chassis Terminal
	Equipotentiality
	On (Supply)
	Off (Supply)
	In Position of a Bi-Stable Push Control
	Out Position of a Bi-Stable Push Control
	Caution: Risk of Electric Shock
	Caution: Hot Surface
	Caution: Risk of Danger
	Warning: Laser Radiation
	Caution: Spinning Blades May Cause Harm

Chapter 2 Safety

Precautions of a general nature should be gathered here. Wherever possible, however, safety warnings, cautions, and notes should only appear immediately before the instructions to which they apply (versus being listed in this section).



DANGER!



The Thorlabs Spectrum Analyzer Controller, SA201, must be powered off, unplugged from the AC input source, and disconnected from any piezo elements prior to replacing the fuse or removal of the cover. Failure to do so may cause **SERIOUS INJURY** to the user, since high voltages exist within the unit.



WARNING!



Complete the setup of the equipment before energizing the laser.



WARNING!



When the laser is energized, always use the proper eye protection relative to the laser used in the setup. When an invisible laser source is used, use extra caution during alignment.

Chapter 3 Overview

Thorlabs' Scanning Fabry-Perot (FP) Interferometers are spectrum analyzers that are ideal for examining fine spectral characteristics of CW lasers. The controller generates a voltage ramp, which is used to scan the separation between the two cavity mirrors. The controller provides adjustment of the ramp voltage and scan time, allowing the user to choose the scan range and speed. An offset control is provided to allow the spectrum displayed on the oscilloscope to be shifted right or left. Another convenient feature of the controller is a zoom capability that provides a 1X, 2X, 5X, 10X, 20X, 50X and 100X increase in the spectral display resolution. The output TTL level trigger allows the user to externally trigger an oscilloscope on either the beginning or midpoint of the ramp waveform.

The SA201 also includes a high precision photodetector amplifier circuit used to monitor the transmission of the cavity. The amplifier provides an adjustable transimpedance gain of 10,000, 100,000, and 1,000,000 V/A when driving a high impedance load, such as an oscilloscope. Using the output sync signal from the controller, an oscilloscope can be used to display the spectrum of the input laser. The detector circuitry incorporates a blanking circuit, which disables the photodiode response during the falling edge of the saw tooth waveform.

3.1. Parts List

Below is a list of all components shipped with the SA201 Spectrum Analyzer Controller.

- SA201 Spectrum Analyzer Controller
- Operating Manual
- 120 VAC US Power Supply Line Cord (with the SA201) or 230 VAC Power Supply Line Cord for Europe (with the SA201-EC)
- 125 mA Fuse for use at 230 VAC operation (250 mA fuse installed in unit)

3.2. Compatible Fabry-Perot Scanning Heads

This product has been designed to be used with one of our SA200, SA210, or SA30 Series Scanning Fabry-Perot Interferometers. Below is a list of available heads. Note this list is subject to change without notice. Please visit the website for the most current information.

Item #	Description
SA200-2B	290 - 335 nm; 520 - 545 nm 1.5 GHz FSR
SA200-3B	350 - 535 nm, 1.5 GHz FSR
SA200-5B	535 - 820 nm, 1.5 GHz FSR
SA200-8B	820 - 1275 nm, 1.5 GHz FSR
SA200-12B	1275 - 2000 nm, 1.5 GHz FSR
SA200-18C	1800 - 2500 nm, 1.5 GHz FSR
SA200-30C	3000 - 4400 nm, 1.5 GHz FSR

Item #	Description
SA210-3B	350 - 535 nm, 10 GHz FSR
SA210-5B	535 - 820 nm, 10 GHz FSR
SA210-8B	820 - 1275 nm, 10 GHz FSR
SA210-12B	1275 - 2000 nm, 10 GHz FSR
SA210-18C	1800 - 2500 nm, 10 GHz FSR
SA30-52	488 - 545 nm, 1.5 GHz FSR

Previous generations of scanning Fabry-Perot interferometers are also compatible with this product. These items are listed below.

Item #	Description
SA200-7A	780 - 930 nm, 1.5 GHz FSR
SA200-9A	900 - 1100 nm, 1.5 GHz FSR
SA200-14A	1450 - 1625 nm, 1.5 GHz FSR
SA200-18B	1800 - 2500 nm, 1.5 GHz FSR
SA200-30B	3000 - 4400 nm, 1.5 GHz FSR

Item #	Description
SA210-5A	525 - 650 nm, 10 GHz FSR
SA210-7A	780 - 930 nm, 10 GHz FSR
SA210-9A	900 - 1100 nm, 10 GHz FSR
SA210-12A	1250 - 1400 nm, 10 GHz FSR
SA210-14A	1450 - 1625 nm, 10 GHz FSR
SA210-18B	1800 - 2500 nm, 10 GHz FSR

Chapter 4 Description

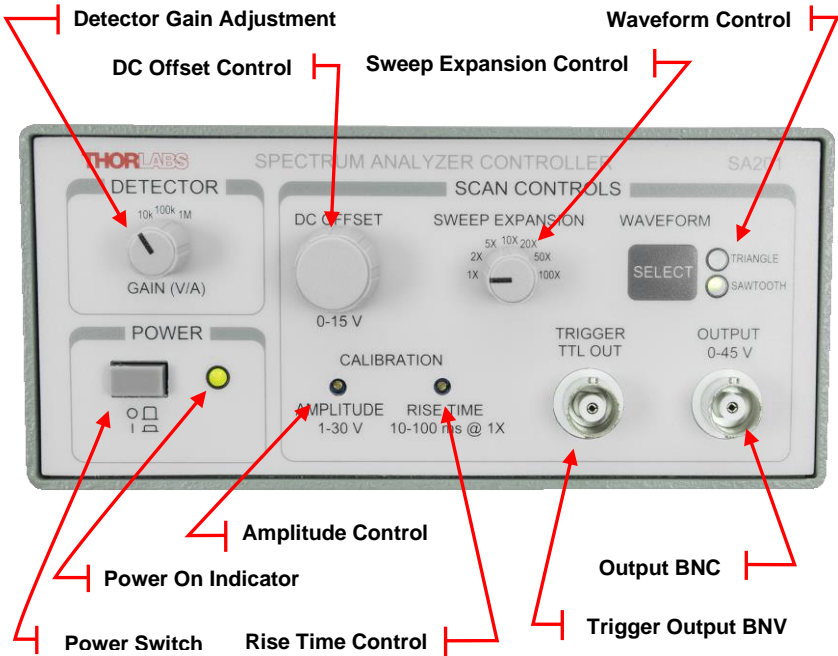


Figure 1 SA201 Front Panel Controls

4.1. Detector Gain Adjustment

The SA201 includes a built in photodiode amplifier circuit. This amplifier is designed specifically to operate with the detector provided with the SA200 series Fabry-Perot Interferometer, allowing the user to monitor the transmission of the cavity. While any photodetector may be connected to the amplifier the specifications, listed in Chapter 3, apply only to detectors supplied with the SA200 series. The amplifier provides a transimpedance gain (current to voltage gain) of 10K, 100K, and 1M V/A while driving a Hi-Z load, such as an oscilloscope. For better noise and performance characteristics it is recommended that a 50 W coax cable with a 50 Ω terminating resistor be used. The photodetector input and output BNC's are located on the rear panel.

4.2. DC Offset Control

The DC Offset provides a continuously adjustable offset voltage over the range of 0 to 15 V using a 10-turn potentiometer. This offset adds directly to the ramp signal. The DC offset control is used to adjust the waveform from left to right across an oscilloscope viewing window, without affecting the calibration of the cavity.

4.3. Sweep Expansion Control

The sweep expansion provides a zoom capability to increase the spectral display resolution by a factor of 1x, 2x, 5x, 10x, 20x, 50x, and 100x. This is achieved by scaling the ramp rise time to the sweep expansion.

4.4. Waveform Control

The SA201 allows the user to select between a saw-tooth and triangular waveform. The saw-tooth waveform is desirable for most applications; however the triangle waveform is useful for cavity alignment. The SA201 will default to the saw-tooth waveform during the system power-up. To change the waveform, simply press the 'WAVEFORM SEL' button. The selected waveform is indicated by the illuminated symbol to the right of the waveform select button.

4.5. Power Switch

The power switch is used to toggle the unit on and off.

4.6. Power On Indicator

The power on LED will light when the unit is powered up.

4.7. Amplitude Control

The amplitude control allows the user to adjust the ramp amplitude from 1 to 30 V peak to peak using a 10-turn trim pot. Note, the ramp signal is added to the DC offset. This means that when the offset is set to 0V, the ramp will start a 0V and increase to the amplitude limit setting. The amplitude is used to determine how far the mirror will be scanned, or to set the spectral range of the optical head.

4.8. Rise Time Control

The rise time control allows the user to continuously adjust the scan rate from 0.01 to 0.1 s using a 10-turn trim pot. Note the rise time setting may be scaled by the sweep expansion setting. For example: If the scan rate is set to 0.05 s and the sweep expansion is adjusted from 1x to 100x then the scan rate will adjust to 5 s. The scaling error is typically less than $\pm 0.5\%$, providing excellent measurement capabilities.

4.9. Trigger Output BNC

This trigger output signal may be used to externally trigger the oscilloscope. The trigger is capable of driving 50 Ω terminated cables, as well as Hi-Z loads such as oscilloscopes. The trigger will provide an edge on the beginning and middle of the scanning ramp. See Figure 2 below.

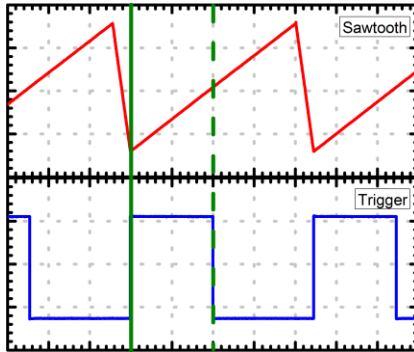


Figure 2 Trigger Logic

4.10. Output BNC

The output BNC is used to drive the SA200 scanning piezos from 1 to 45 V. The output is capable of driving 0.6 μF piezo loads at a ramp rate of 1 ms over the full voltage range. The output current is internally limited to prevent damage to the output drive. Note: the output performance specifications assume a Thorlabs Fabry-Perot Interferometer module is connected.

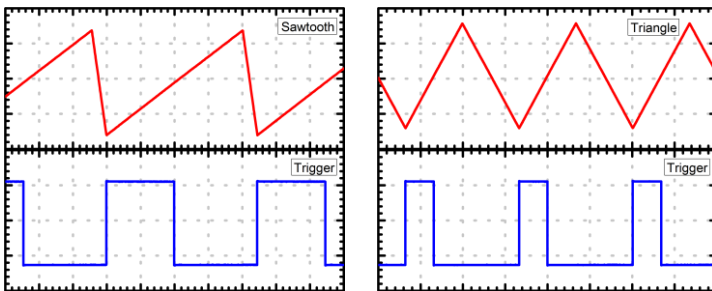


Figure 3 Saw-tooth Waveform (left), Triangle Waveform (right)

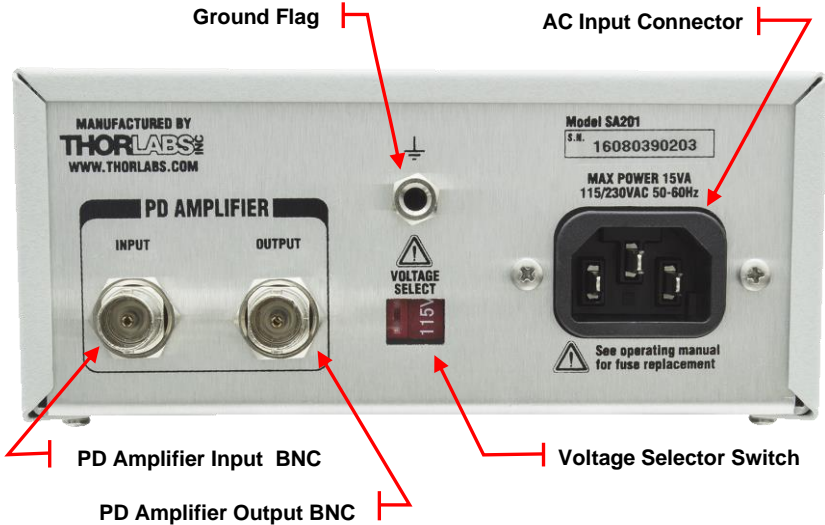


Figure 4 SA201 Rear Panel Connections

4.11. Ground Plug

This ground plug is for use as a general-purpose ground power connection. It is connected directly to the earth ground connection of the input power plug.

4.12. AC Input Connector

This is the line voltage input connection. **IMPORTANT:** The unit is configured for 100/115 VAC, 50 – 60 Hz from the factory. To operate at 230 VAC see Section 5.3.

4.13. PD Amplifier Input BNC

This input BNC is used to interface the photodetector, provided with the SA200 scanning heads, to the amplifier circuit. The photodiode amplifier is configured to operate with the Thorlabs supplied photodetectors; however it is possible to operate user supplied photodetectors. To do so, the BNC center contact must be connected to the photodetector cathode and the BNC shell must be connected to the photodiode anode (unbiased operation). If a biased detector is to be used the BNC shell must be connected to the bias ground and the bias voltage must be negative for the circuit to operate properly.

**NOTE**

The PDAVJ5 that is recommended for use along with the SA200-30C should not be connected to this input. Connecting to the amplifier will cause damage to the instrument.

4.14. PD Amplifier Output BNC

This BNC is the amplifier output and may be connected directly to an oscilloscope to view the cavity spectrum. The amplifier gain will be set using the front panel 'DETECTOR' control knob. The amplifier output includes a 50 Ω series resistor to minimize noise when operating with a 50 Ω coax cable. For best results, a 50 Ω load resistor is recommended at the oscilloscope. Note, the amplifier gain will be halved with a 50 Ω load connected.

4.15. Voltage Selector Switch

The voltage selector switch allows the user to select the input line voltage. The factory default setting is 100/115 VAC as shown in Figure 4. To operate at 230 VAC, this switch will have to be moved to the 230 V position. The line fuse will also need to be changed to properly protect the unit. See sections 5.2 below for detailed instructions.

Chapter 5 Operation

5.1. PD Blanking Circuit

The detector amplifier includes a blanking circuit, which blocks any photodetector response during the falling edge of the saw tooth waveform. This is very useful when triggering on the photodiode spectral response, because unwanted signals while the cavity resets will be removed. The blanking is not available when using the triangle waveform, since it is useful to see the rising and falling response overlapped during system alignment. This feature may be disabled as described below:



IMPORTANT



**Before continuing, disconnect the scanning head or any piezo device from the SA201 output. Disconnect the power cord.
Do not open the unit if the power cord is connected.**

1. Remove the two screws securing the enclosure cover with a Phillips head screwdriver. The screws are located on the bottom side, rear corners of the unit. Do not lose the screws.
2. Carefully remove the cover by sliding toward the rear of the unit.
3. Locate the JP3 header. It is positioned in front of the heat sink and will have a shorting jumper on pin 1.

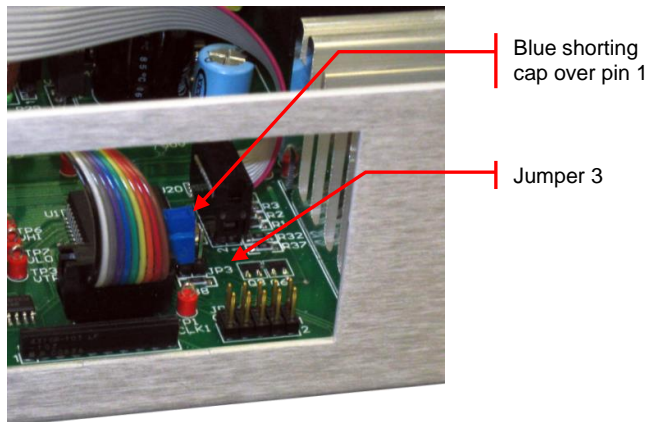


Figure 5 Shorting Cap and Jumper 3

4. Remove the shorting jumper and place across (shorting) the JP3 pins to disable the blanking circuit. The default setting will be blanking enabled. The jumper will not be shorting the pins.
5. Replace the enclosure cover and secure with the enclosure screws.

5.2. Replacing the Fuse



DANGER!



The Thorlabs SA201 Spectrum Analyzer Controller must be powered off, unplugged from the AC input source, and disconnected from any piezo elements prior to replacing the fuse or removal of the cover. Failure to do so may cause **SERIOUS INJURY** to the user, since high voltages exist within the unit.

5.2.1. Materials Needed

- SA201 Operating Manual – The most recent version of this operating manual will be available on the Thorlabs web site.
- 250 mA Type 'T' Slow Blow Fuse – The 250 mA fuse is installed from the factory. This must be installed when operating the unit at 100/115 VAC.
- 125 mA Type 'T' Slow Blow Fuse – The 125 mA fuse is required for 230 V operation only. Thorlabs supplies a 125 mA fuse with all of its SA201 units and must be installed when operating at 230 VAC.
- Phillips Head Screwdriver (#2 Preferred) – We do not recommend using electrically powered screwdrivers.

5.2.2. Fuse Replacement



IMPORTANT



**Disconnect the scanning head or any piezo device from the SA201 output.
Disconnect the power cord.
Do not open the unit if the power cord is connected.**

1. Remove the two screws securing the enclosure cover with a Phillips head screwdriver. The screws are located on the bottom side, rear corners of the unit. Do not lose the screws.
2. Carefully remove the cover by sliding toward the rear of the unit.

3. Locate the fuse box between the input line voltage connector and the transformer.

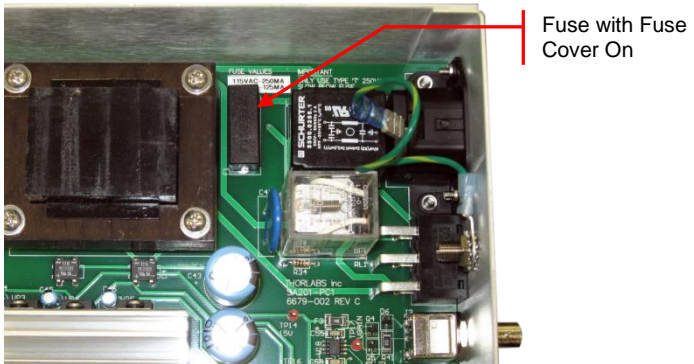


Figure 6 SA201 Fuse with Fuse Cover

4. Remove the fuse cover and slide the old fuse out.
5. Install the new fuse into the fuse cover and place back into the fuse box. (250 mA for 100/115 VAC and 125 mA for 230 VAC)
6. Replace the enclosure cover and secure with the enclosure screws.

5.3. Selecting the Line Voltage

1. Replace the line fuse as described above.
2. Locate the voltage selector switch on the rear panel. See Figure 4 above.
3. Switch to the appropriate line voltage.
4. Install the appropriate line cord.

5.4. Cleaning

The SA201 should only be cleaned with a soft cloth and a mild soap detergent or isopropyl alcohol. Do not use a solvent-based cleaner.

Chapter 6 Recommended Setup

6.1. Recommended Setup for SA200 Series Fabry-Perot Interferometers (Except SA200-30C)

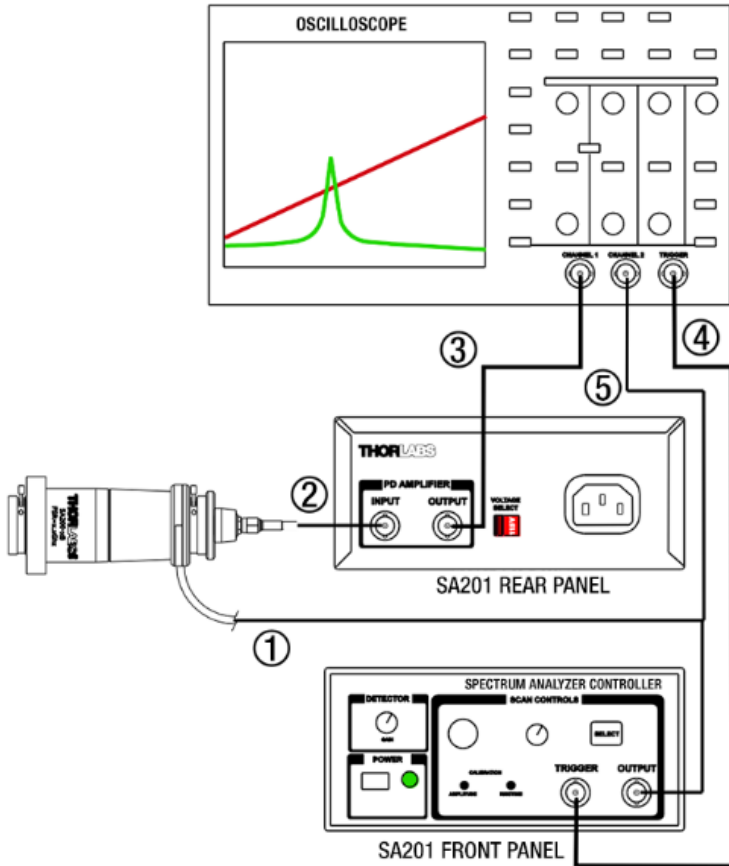


Figure 7 Recommended Setup Diagram (Except SA200-30C)

6.2. Recommended Setup for SA200-30C Fabry-Perot Interferometer

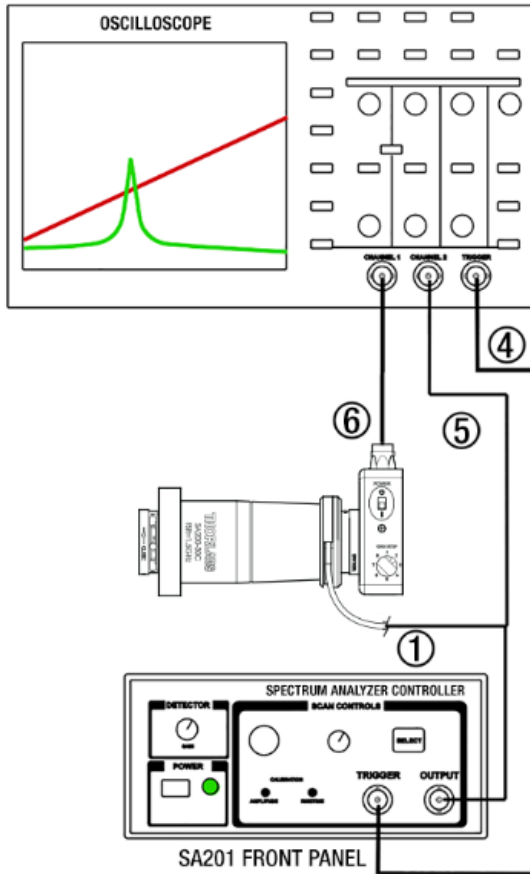


Figure 8 Recommended Setup Diagram for SA200-30C

Connection	Description
1	Controller (BNC) to Piezo (Cable is Attached to FP Interferometer)
2 ¹	Photodiode (SMA) to Controller (BNC) (Included with FP Interferometer)
3 ¹	Amplified Photodiode Output (BNC) to Oscilloscope (Not Included)
4	Trigger Output of Controller (BNC) to Oscilloscope (Not Included)
5	Optional Connection that Allows the User to Monitor the Signal used to Drive the Piezoelectric Transducers (Not Included)
6 ²	PDAVJ5 Output (BNC) to Oscilloscope (Detector and Cable Not Included)

¹ This connection is not part of the setup for the SA200-30C.

² This connection is part of the setup only for the SA200-30C.

Chapter 7 Specifications

7.1. Output Characteristics

Item #	SA201
Waveforms	Sawtooth / Triangle
Default Waveform	Saw tooth
Saw tooth Fall Time	1 ms Typical
Output Voltage Range	1 to 45 V (Offset + Amplitude)
Max Supply Current ³	15 mA
Short Circuit Current ⁴	26 mA Max
Short Circuit Duration ⁴	Continuous
Offset Adj. Range	0 to 15 VDC
Amplitude Adj. Range	1 to 30 V
Rise Time Adj. Range ⁵	0.01 to 0.1 s @ 1X Sweep Exp. 1 to 10 s @ 100X Sweep Exp.
Sweep Expansion Settings	1X, 2X, 5X, 10X, 20X, 50X, 100X
Sweep Scale Error ⁶	±0.5%
Output Noise ⁷	1 mV _{RMS} (~6.6 mV _{PP})

7.2. Trigger Characteristics

Item #	SA201
Trigger Output Voltage	TTL levels
VOH (R _L = 50 W)	2 V Min
VOL (R _L = 50 W)	0.5 V Max
Trigger Load Impedance	50 W / Hi-Z
Trigger Rising Edge ⁸	Ramp Start
Trigger Falling Edge ⁸	Ramp Midpoint

³ Achieved during the saw tooth waveform fall time. This is calculated by

$$I \text{ (mA)} = \frac{(C_{piezo}(\mu\text{F}))(\Delta V_{max})}{\Delta t_{fall}}$$

⁴ The output drive amplifier will current limit the load to 26 mA max. Although the unit may operate continuously under these conditions, it is not recommended since the unit will heat up causing stress to the electronics.

⁵ The rise time adjustment range for each sweep setting is as follows:
=(0.01 x "Sweep Expansion Setting")" to "(0.1x"Sweep Expansion Setting")

⁶ Defined as the scaling error between 1X and any other gain settings (ex. 2X ± 0.5%).

⁷ Measured with SA200 series scanning head connected to output.

⁸ 'Ramp' refers to the rising, or scanning, edge of the 'Output' waveform.

7.3. Photodiode Amplifier Characteristics⁹

Item #	SA201
Gain Steps	0, 10, 20 dB
Transimpedance Gain (Hi-Z)	10K, 100K, 1M V/A
Transimpedance Gain (50 Ω)	5K, 50K, 500K V/A
Gain Error ¹⁰	$\pm 0.1\%$ @ 10K ($\pm 0.12\%$) $\pm 0.12\%$ @ 100K ($\pm 0.15\%$) $\pm 0.14\%$ @ 1M ($\pm 0.3\%$)
Output Impedance	50 Ω
Load Impedance	50 Ω /Hi-Z
Output Voltage (Hi-Z load)	0 - 10 V Min
Output Voltage (50 Ω load)	0 - 5 V Min
Max Output Current ¹¹	100 mA
Bandwidth ¹¹	250 KHz
Noise (RMS) ¹¹	<0.1 mV @ 10K 0.2 mV @ 100K 1.5 mV @ 1M
Offset ¹¹	± 1 mV @ 10K ± 5 mV @ 100K ± 20 mV @ 1M

⁹ These characteristics are not for the SA200-30C and the recommended PDAVJ5 detector.

¹⁰ The gain error does not apply when using a 50 Ω load since the user-installed output terminator will probably have a resistance tolerance greater than the gain errors above. Also note that the 50 W output series resistance is 49.9 W $\pm 1\%$. This will also factor into gain error when using a 50 Ω load.

¹¹ Test performed with a 50 Ω terminator and a 6' (~1.8 m) 50 Ω coax cable.

7.4. Physical Features

Item #	SA201
Dimensions (W x H x D)	5.8" x 2.8" x 12.5" (147 mm x 71 mm x 317.5 mm)
Input and Output Connectors	BNCs
Offset Control	10-Turn Potentiometer
Amplitude Control	10-Turn Trim pot
Rise Time Control	10-Turn Trim pot
Sweep Expansion Control	7-Position Rotary Switch
Photodiode Gain Control	3-Position Rotary Switch
Waveform Select	Pushbutton w/ Illuminated Indicators
PD Amplifier Features	Blanking with Sawtooth Waveform Falling Edge
Operating Temperature	10 to 40 °C
Storage Temperature	0 to 85 °C

7.5. Power Supply

Item #	SA201
Supply Type	Linear
Voltage Selection	Switch Selectable between 115 / 230 VAC
Input Voltage	100 / 115 / 230 VAC
Line Frequency	50 - 60 Hz
Input Power	15 W Max
Fuse Ratings	250 mA @ 100 / 115 VAC, 125 mA @ 230 VAC
Fuse Type	Slow Blow Type 'T'

Chapter 8 Regulatory

As required by the WEEE (Waste Electrical and Electronic Equipment Directive) of the European Community and the corresponding national laws, Thorlabs offers all end users in the EC the possibility to return “end of life” units without incurring disposal charges.

- This offer is valid for Thorlabs electrical and electronic equipment:
- Sold after August 13, 2005
- Marked correspondingly with the crossed out “wheelie bin” logo (see right)
- Sold to a company or institute within the EC
- Currently owned by a company or institute within the EC
- Still complete, not disassembled and not contaminated



As the WEEE directive applies to self contained operational electrical and electronic products, this end of life take back service does not refer to other Thorlabs products, such as:

- Pure OEM products, that means assemblies to be built into a unit by the user (e.g. OEM laser driver cards)
- Components
- Mechanics and optics
- Left over parts of units disassembled by the user (PCB's, housings etc.).

If you wish to return a Thorlabs unit for waste recovery, please contact Thorlabs or your nearest dealer for further information.

Waste Treatment is Your Own Responsibility


If you do not return an “end of life” unit to Thorlabs, you must hand it to a company specialized in waste recovery. Do not dispose of the unit in a litter bin or at a public waste disposal site.

Ecological Background

It is well known that WEEE pollutes the environment by releasing toxic products during decomposition. The aim of the European RoHS directive is to reduce the content of toxic substances in electronic products in the future.

The intent of the WEEE directive is to enforce the recycling of WEEE. A controlled recycling of end of life products will thereby avoid negative impacts on the environment.

Chapter 9 Declaration of Conformity

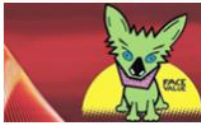

THORLABS
www.thorlabs.com

EU Declaration of Conformity

in accordance with EN ISO 17050-1:2010

We: Thorlabs Inc.

Of: 56 Sparta Avenue, Newton, New Jersey, 07860, USA

in accordance with the following Directive(s):

2014/35/EU	Low Voltage Directive (LVD)
2014/30/EU	Electromagnetic Compatibility (EMC) Directive
2011/65/EU	Restriction of Use of Certain Hazardous Substances (RoHS)

hereby declare that:

Model: **SA201 & SA201-EC**

Equipment: **Controller for Scanning Fabry-Perot Interferometers**

is in conformity with the applicable requirements of the following documents:

EN 61010-1	Safety Requirements for Electrical Equipment for Measurement, Control and Laboratory Use.	2010
EN 61326-1	Electrical Equipment for Measurement, Control and Laboratory Use - EMC Requirements	2013

and which, issued under the sole responsibility of Thorlabs, is in conformity with Directive 2011/65/EU of the European Parliament and of the Council of 8th June 2011 on the restriction of the use of certain hazardous substances in electrical and electronic equipment, for the reason stated below:

does not contain substances in excess of the maximum concentration values tolerated by weight in homogenous materials as listed in Annex II of the Directive

I hereby declare that the equipment named has been designed to comply with the relevant sections of the above referenced specifications, and complies with all applicable Essential Requirements of the Directives.

Signed:  On: 09 April 2018

Name: Ann Strachan

Position: Compliance Manager

EDC - SA201 & SA201-EC -2018-04-09



Chapter 10 Thorlabs Worldwide Contacts

For technical support or sales inquiries, please visit us at www.thorlabs.com/contact for our most up-to-date contact information.



USA, Canada, and South America

Thorlabs, Inc.
sales@thorlabs.com
techsupport@thorlabs.com

Europe

Thorlabs GmbH
europe@thorlabs.com

France

Thorlabs SAS
sales.fr@thorlabs.com

Japan

Thorlabs Japan, Inc.
sales@thorlabs.jp

UK and Ireland

Thorlabs Ltd.
sales.uk@thorlabs.com
techsupport.uk@thorlabs.com

Scandinavia

Thorlabs Sweden AB
scandinavia@thorlabs.com

Brazil

Thorlabs Vendas de Fotônicos Ltda.
brasil@thorlabs.com

China

Thorlabs China
chinasales@thorlabs.com



THORLABS

www.thorlabs.com
