

## INF0091T – PSY35x ScienceDesk Side Shelf - Assembly Instructions

This information sheet details the assembly instructions for the PSY35x Side Shelf.

For the continuing safety of the operators of the equipment and for the protection of the equipment itself, the operator should take note of the **Warnings, Cautions and Notes** throughout this leaflet, and, where applicable, in any associated literature.

### 1.1 Fitting the Mounting Brackets

Referring to Fig. 1.1:

- 1) Position the brackets on the underside of the shelf as shown below, such that the side face is outermost.
- 2) Using the six posi head screws supplied, attach the brackets to the shelf.
- 3) Tighten the screws.

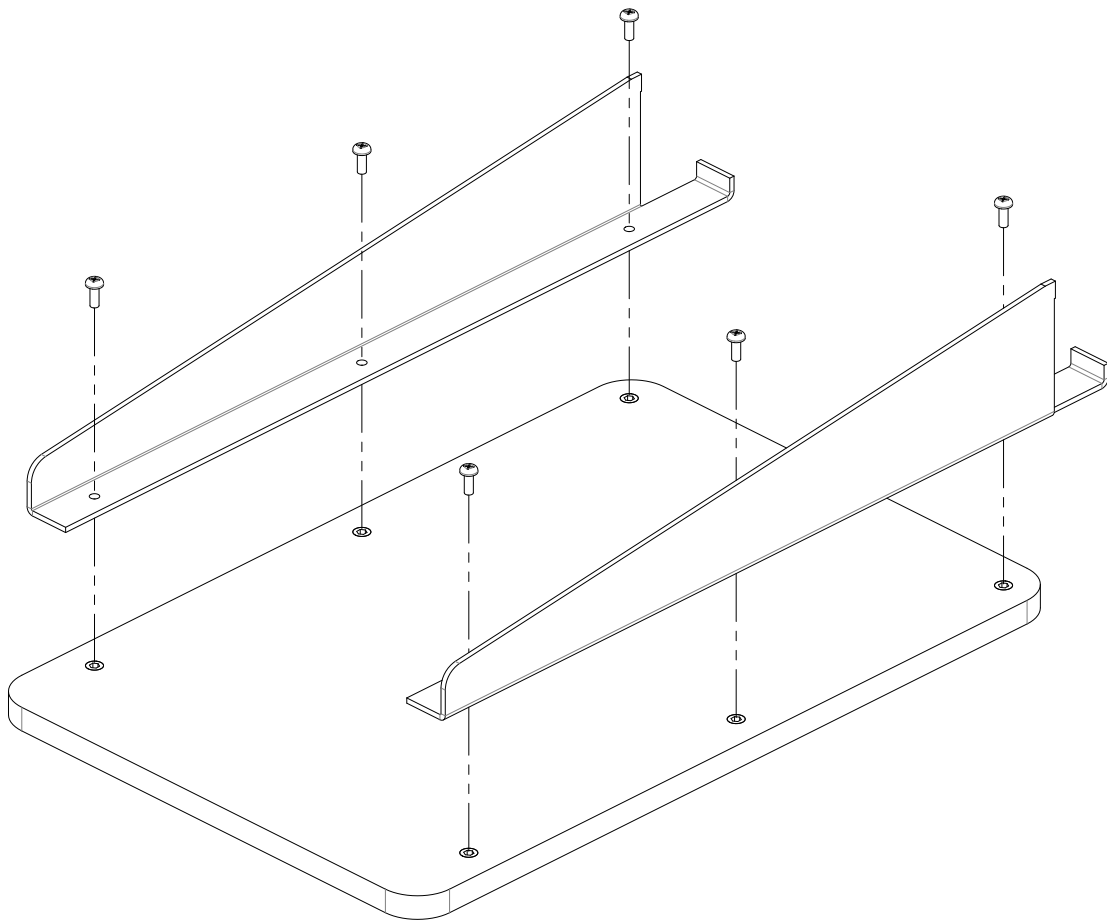


Fig. 1.1 Fitting the attachment brackets

## 1.2 Fitting the Side Shelf

The side shelf is to be fitted to the upper side rails of the desk and is designed to support a mass up to 50kg.



Referring to Fig. 1.2:

- 1) Position the shelf on the upper rail of the desk.
- 2) Locate the shelf clips over the rail and position for user convenience.
- 3) Ensure no wires are trapped under the shelf or the mounting brackets.



Fig. 1.2 Fitting the side shelf

